

Specification

Incuwater-Clean™

A5219

HS-No.:	38220000
Storage:	2-8°C
	<ul style="list-style-type: none"> • Disinfectant solution for CO₂-incubator water baths. • for the prevention of microbial growth in incubator water baths. • 100X concentrated solution. Use 50 ml per 5 liter of incubator water bath.
	<ul style="list-style-type: none"> • Incuwater-Clean™ is a registered trademark of AppliChem GmbH.
<p>Comment</p> <p>The water required to create the humidity is a source of contamination which disperses in the incubator. In order to disinfect the water we recommend Incuwater-Clean™ solution, which contains a disinfectant that does not cause damage to the stainless steel tray, is non-toxic, non-volatile, and extremely effective. The water should be replaced with sterile water every two to four weeks, adding 50 ml of Incuwater-Clean™ per 5 liters of water. This product can be applied in normal water baths too.</p> 	